


Procedure for connecting M5 cables to connectors.


The M5 cable supplied with our photo-detector system has male pins on both ends. Therefore the connection procedure is the same for the FDA (amplifier) and FDH (head).

doriclenses.com


1 The connectors have a key.

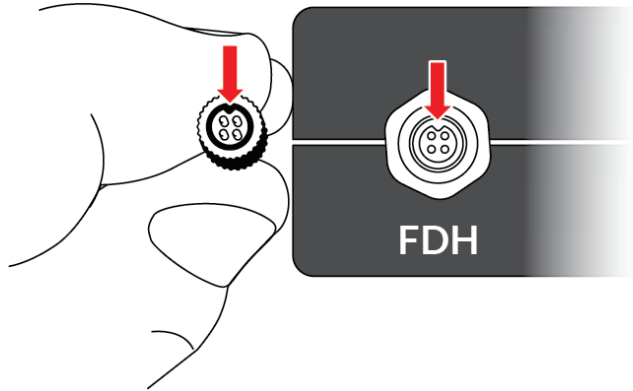


3 Without rotating the cable, continue inserting it into connector until it reaches a stop. If the keys are well aligned, the threads should be in part inside the connector.

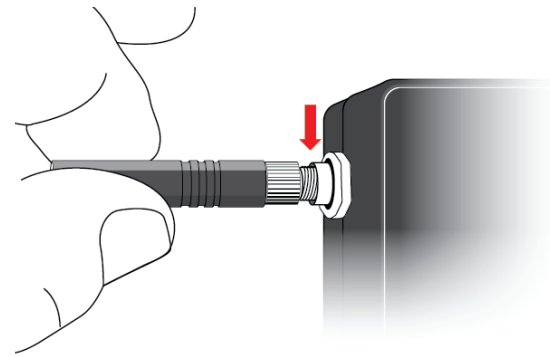

2 Make sure that the keys on both ends are aligned and gently insert the cable.

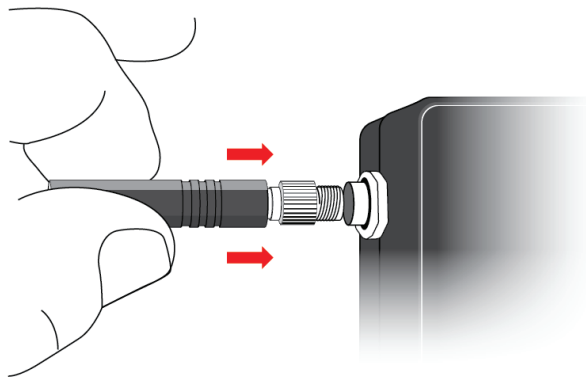

4 Turn clockwise the metal barrel of cable connector to tighten it.

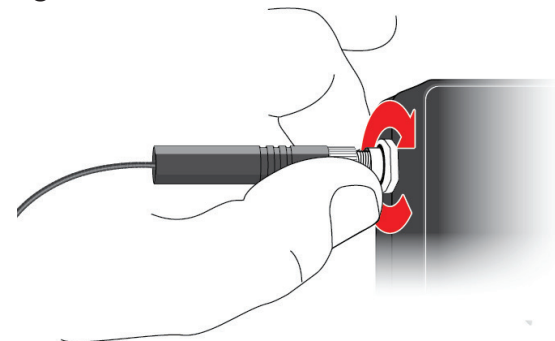

Note: While tightening the connector, make sure to turn only the metal barrel, NOT the whole cable.

doric

5 To disconnect the cable, please follow the same procedure in reversed order.